



## LEAVES NO VISIBLE RESIDUE REDUCING THE NEED FOR FREQUENT RINSING



### Klercide™ Cleanroom Quat\* Blended with WFI

A quaternary ammonium compound solution specifically designed for Pharmaceutical cleanrooms

**A ready-to-use sanitizer ideal for routine use on hard surfaces. The product is efficacious at a low concentration, ensuring the product leaves minimal residues. This makes it ideal for high grade cleanrooms and critical manufacturing areas.**

Klercide Cleanroom Quat is available in multiple formats, including a one gallon, capped, ready-to-use product and a 32 oz. ready-to-use product.

Klercide Cleanroom Quat is:

- ▲ A quaternary ammonium compound with broad spectrum efficacy
- ▲ Available double or triple bagged
- ▲ Supplied in our proprietary and validated SteriShield Delivery System

The SteriShield Delivery System (SDS) features:

- ▲ A closed trigger system providing a completely sealed system
- ▲ Protection for the integrity of the contents throughout use
- ▲ Validated in-use shelf life of six months
- ▲ Dispensing of the full volume of product
- ▲ A fully adjustable spray pattern to ensure optimum surface coverage

#### PRODUCT QUALITY ASSURANCE

Manufactured in our state-of-the-art facilities to the principles of GMP, Klercide Cleanroom Quat is safe for use even in critical aseptic core manufacturing areas.



\*Same product available under a different name (Klercide Low Residue Quat) outside of US and Canada.

# Klercide™ Cleanroom Quat\*

Blended with WFI



A quaternary ammonium compound solution specifically designed for Pharmaceutical cleanrooms

## Why use Klercide Cleanroom Quat for routine sanitization?

### OPERATOR SAFETY

Klercide Cleanroom Quat is effective at a lower concentration, which reduces the likelihood of sticky floors, slips and falls as well as unfavorable cross-reactions which can cause hazardous vapors.

### REDUCED COST OF USE

Klercide Cleanroom Quat eliminates the need for frequent or daily residue management which may help increase production time and reduce costs related to residue management such as labor and time associated with additional cleaning.

### COMPATIBLE

Klercide Cleanroom Quat can be used on all types of hard surfaces in cleanrooms and controlled environments, isolators, equipment and production vessel exteriors. It is compatible with other Ecolab products including Klercide Low Residue Peroxide - and is therefore suitable for rotational use.

### DESIGNED FOR CLEANROOMS

Klercide Cleanroom Quat is manufactured in Ecolab's ISO 9001 and ISO 14001 accredited facilities to the principals of GMP. These manufacturing conditions ensure Klercide Cleanroom Quat is manufactured to the highest quality and is compliant for use in your critical areas.

Klercide Cleanroom Quat is:

- ▲ Made with a quaternary ammonium compound
- ▲ Blended with WFI
- ▲ Endotoxin tested to an approved pharmacopoeia method
- ▲ Packed and double or triple bagged in ISO Class 5 cleanroom
- ▲ 0.2 micron filtered and aseptically filled under unidirectional flow
- ▲ Supported with Certificates of Analysis and Sterility for every batch
- ▲ 1-year shelf life
- ▲ Supported with full documentation

## MICROBIOLOGICAL TEST RESULTS

EFFICACY	CONTACT TIME
Non-Food Contact Sanitizer	5 min

Please see label for specific claim

### WORLDWIDE HEADQUARTERS

1 Ecolab Place  
St. Paul, MN 55102  
USA

[www.ecolab.com/lifesciences](http://www.ecolab.com/lifesciences)  
[www.ecolab.com/pharmaceutical](http://www.ecolab.com/pharmaceutical)

© 2019 Ecolab USA Inc. All rights reserved. 20FEB19/NA

PHARMACEUTICAL

## Quality. Compliance. Safety.

### APPLICATION AREA

For use on floors, walls and hard surfaces.

### INSTRUCTIONS FOR USE

**Spray:** For optimum results hold bottle upright and spray from a distance of 6" to 8". Always close the nozzle after use.

**Wetting:** Apply solution with mop, disposable wiper, coarse spray device or by soaking so as to wet all surfaces thoroughly. Allow surfaces to remain wet for a minimum of 5 minutes and then either allow to air dry or if desired wipe dry with a (sterilized) cloth or wiper. See label for complete directions for use.

### INGREDIENTS

Didecyl dimethyl ammonium chloride and Water For Injection.

### REGISTRATION DETAILS

EPA Reg. No. 1677-257

### PACKAGING INFORMATION

Cleanroom Quat Blended with WFI

UNIT SIZE	UNIT OF SALE
32 oz spray	6 x 32 oz
1 Gallon capped	4 X 1 gal.

FOR PRODUCT ORDER CODES AND ADDITIONAL INFORMATION, PLEASE CONTACT YOUR ACCOUNT MANAGER OR VISIT OUR WEBSITE

### SHELF LIFE

Unopened, this product has a one year shelf life.

ALWAYS READ THE LABEL AND PRODUCT INFORMATION BEFORE USE.

**ECOLAB®**